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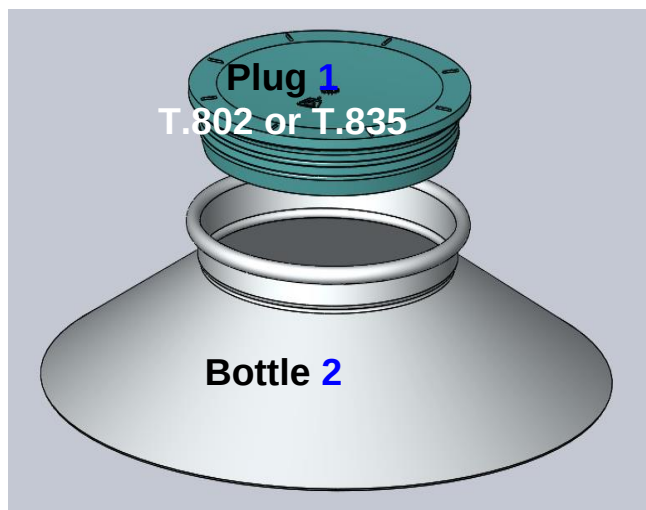
PLUG INSERTION / REMOVAL TOOL

T.802 (item 7319) or T.835 (item 7327)

IMPORTANT: This insertion/removal tool is preset and tested in our premises before despatch. Any major adjustment or dismantling should be made at TOURNAIRE. In case of any doubt, do not hesitate to contact us or send us the tool for adjustment. During the 2 year warranty period, TOURNAIRE can ensure the repair or adjustment of the tool, provided it has not been submitted to abnormal mechanical or chemical constraints nor used improperly. We recommend for instance not to sterilize/ depyrogenize with oxidant substances (ethylene oxide, hydrogen peroxide,..) and rather use autoclave or Dry Heat Sterilization for this purpose.

Closure System T.802 or T.835

Picture A

1 - PREPARATION:

1-1 Check the insertion/removal tool before use:

Before use (picture A), make sure that the insertion/removal tool 3 alone operates properly by pushing downwards the levers: you should not feel any rubbing, unusual resistance or abnormal noise.

In case of malfunction, refer to §4 or contact **TOURNAIRE** for assistance.

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1-2 Preparation of bottle and closure:

- Position the container **2** on a flat surface at a convenient height for the operator (neck of bottle at hips height).
- Place the container close to the perforated stainless steel base **4** (item 7325) - see picture **B**

2 – INSERTION OF THE PLUG:


- After cleaning and eventually sterilizing by autoclave , place the plug **1** on the base **4** for this purpose (picture **C**)
- Place the insert/removal tool over the plug (picture **D**). The tool must be perfectly horizontal centered and in full contact all around the plug (picture **E**).
- Simultaneously push the levers downwards (picture **F**) until you reach the horizontal blocking position and levers are stopped
- Lift the tool and plug keeping levers horizontal and position the tool and plug centered inside the neck of the bottle (picture **G**)
- Press firmly downwards on the tool to insert fully the plug (picture **H**)
- Lift the levers vertically to release the plug (picture **I**).
- Remove the tool and check the conformity of insertion:

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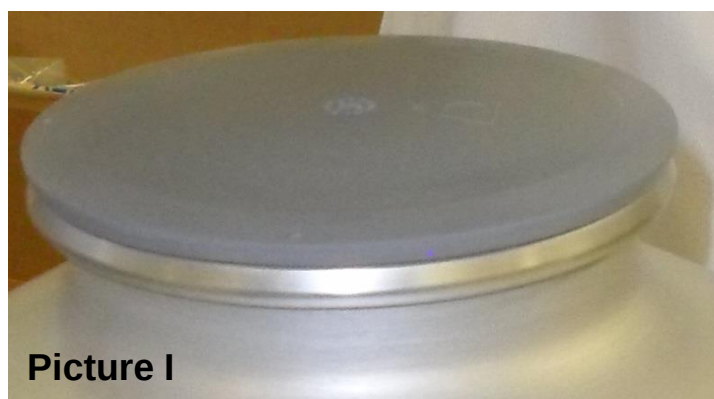the plug **1** must be in full contact all around the neck of the container. (Picture I)

3 - REMOVAL OF THE CAP:

- Repeat every insertion instruction of §2 in reverse order.

4 - TROUBLE SHOOTING:

In any case, first ensure that the plug and the Insert/removal tool are adequate and well prepared – see § 1-1 and 1-2.

1/ too much effort is required:

- Check that there are no foreign bodies, particles or dirt in the tool. If necessary, clean it with a cloth soaked in alcohol.
- Check if the instructions of section 1.2 have been met: proper placement of the cap.

2/ plug does not adhere to the tool:

- Check that the plug is in contact with the bottom of the inner part of the tool.

In case of doubt, do not hesitate to contact Tournaire for assistance

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En cas de mauvais fonctionnement, voir §4 ou contacter **TOURNAIRE S.A.**