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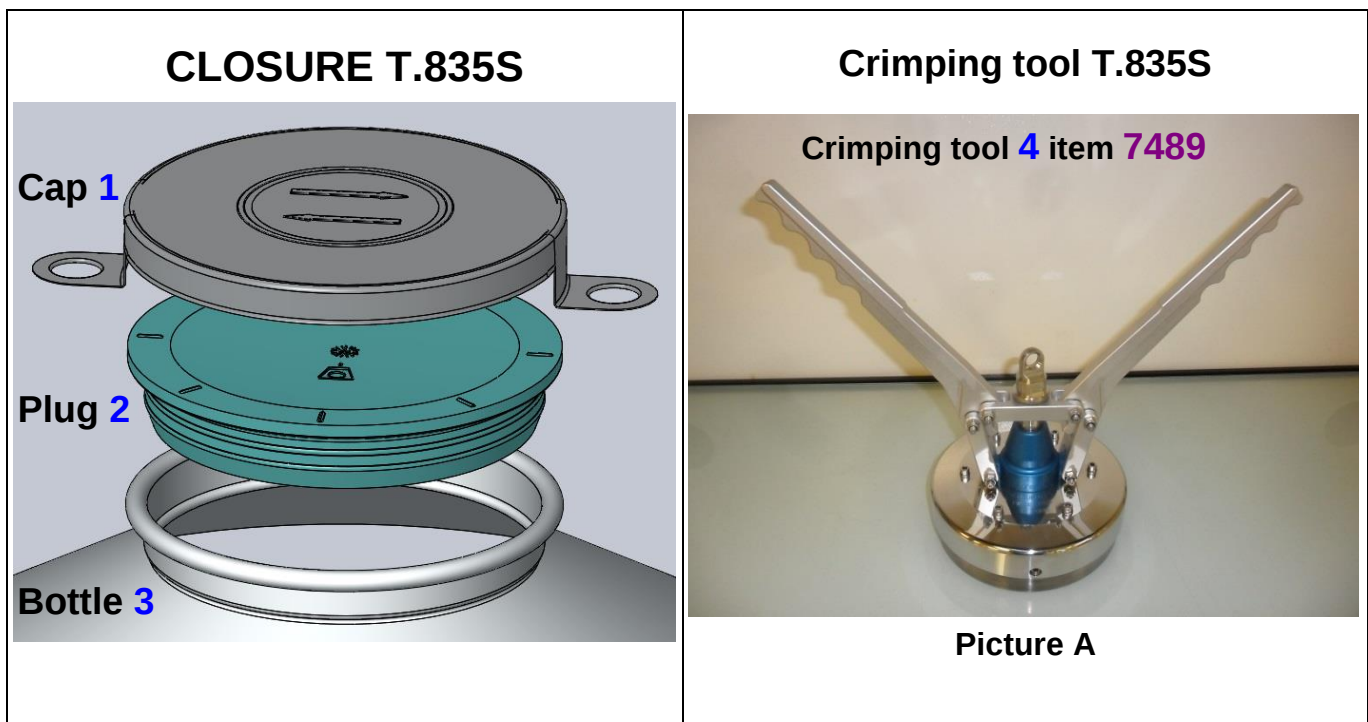
USER'S GUIDE

Sealing / Crimping tool T 835S

IMPORTANT: This crimping tool is pre-set and tested in our premises before despatch. Any major adjustment or dismantling should be made at TOURNAIRE. In case of any doubt, do not hesitate to contact us or send us the tool for adjustment. During the 2 year warranty period, TOURNAIRE can ensure the repair or adjustment of the tool, provided it has not been submitted to abnormal mechanical or chemical constraints nor used improperly. We recommend for instance not to sterilize/ depyrogenize with oxidant substances (ethylene oxide, hydrogen peroxide,...) and rather use autoclave or Dry Heat Sterilization for this purpose.

1 - PREPARATION:

1.1 Check the crimping tool before use :

Before crimping, make sure that the crimping tool 4 alone operates properly by pushing downwards the levers: you should not feel any rubbing, unusual resistance nor abnormal noise. Check also internal surface of the tool and rotating parts: they should be free from foreign objects that could impair the crimping.


In case of malfunction, please check that proper maintenance has been carried out – see **MAINTENANCE FORM NOTI-0028** or contact **TOURNAIRE S.A.** for assistance.

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
1.2 – Preparation of bottle and closure accessories before crimping

Precaution: Avoid direct contact with accessories and protect yourself with suitable gloves.

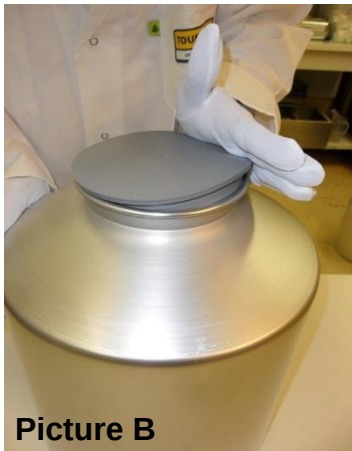- Place the container on a flat surface at a convenient height for the operator (neck of bottle ideally at hips height).
- Place the plug **2** either manually – see picture **B** and **C**, or use the inserting/removal tool **5** - Ref. TOURNAIRE **5314** or **7327** see picture **D**. Make sure that the edge of the plug **2** is in full contact all around the container rolled-on neck.
- Place the aluminium cap **1** on the plug **2** on the top of it (pictures **E**, **F**, **G**), push firmly on the top, make sure that the top of cap covers entirely the plug and is in contact with the plug everywhere.

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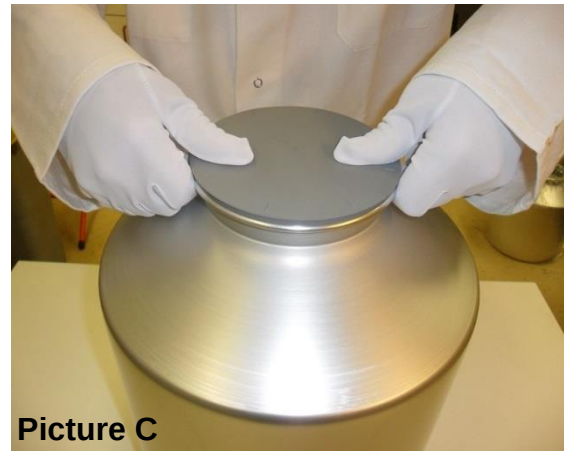
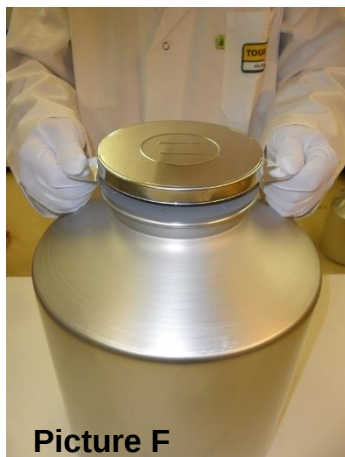
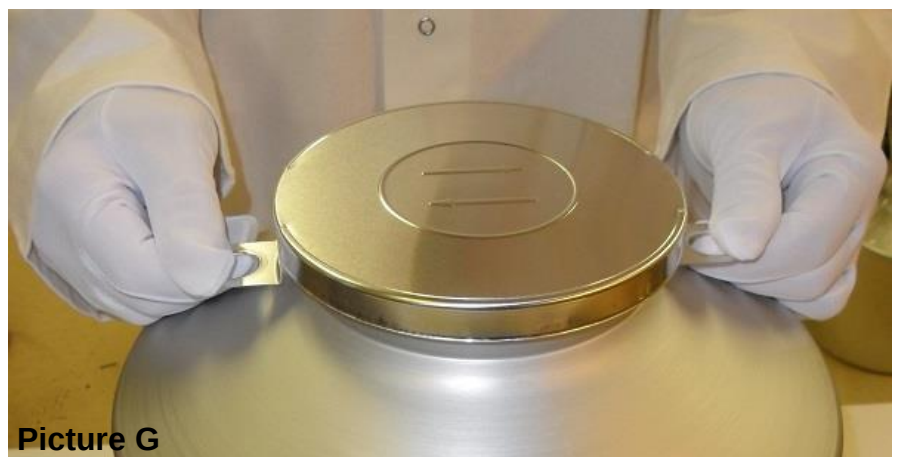
2 - CRIMPING OPERATION:

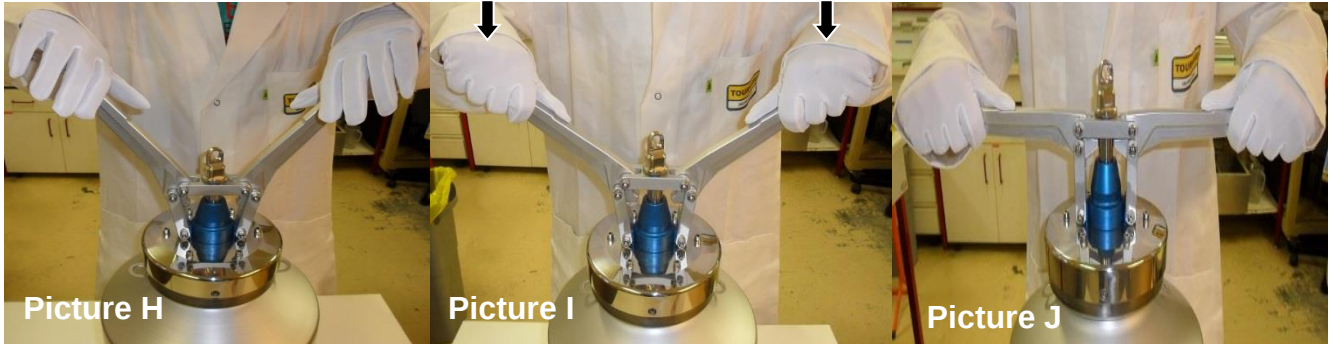- Position the crimping tool on the assembled closure and bottle. Press firmly downwards so that the tool is strictly horizontal and fully in contact with the aluminium cap (picture **H**).
- Simultaneously push the levers downwards (picture **I**) until you reach the horizontal blocking position and levers are stopped (picture **J**). The cap is crimped.
- Lift the levers.
- Remove the crimping tool.

3 - QUALITY CONTROL of the CRIMPING:

The aluminium cap **1** must be entirely rolled back on the container's neck all around the neck circumference, some little folds on the edge of the cap don't reduce the performance of the sealing (picture **K**). It is normal to see little folds on the edge of the cap.

4 - TROUBLE SHOOTING:

In case a defective crimping is noted, first check that plug, cap and tool are properly adjusted – see § 1-1 and 1-2. If cap is blocked or deformed, take it off and use a new one then retry.

1/ too much effort is required for crimping:

- Check the crimping tool is free from any foreign object or dirt. Clean with a tissue or cloth, dipped into alcohol for instance and remove the dirt /object.
- Check that all preparation steps (§1.2) have been thoroughly respected, particularly the full horizontal insertion of the plug and cap, everywhere in contact.
- Repeat crimping operations (see - §2 -) with a new aluminium cap **1**
- Check that the crimping tool has been well maintained (see especially NOTI-0028 §4)

2/ crimping still not conform (Cf. § 3):



In case of malfunction of the crimping tool , doubt or incorrect crimping, please check that proper **MAINTENANCE FORM-NOTI-0028** has been carried out

In case of doubt, do not hesitate to contact TOURNAIRE .

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5 –CAP OPENING INSTRUCTION:

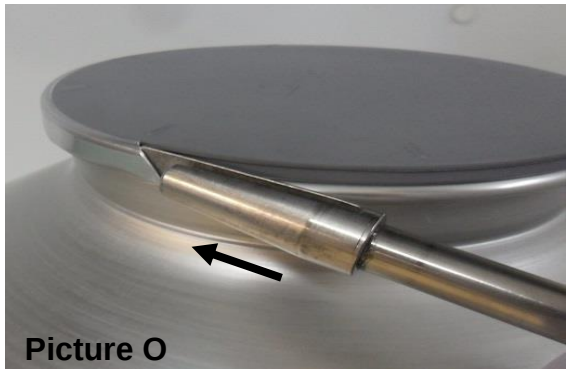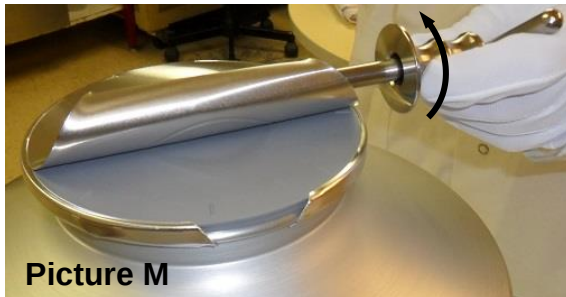- If not easily reachable, lift one of the cap ears upward with the edge of the cap opener **6** and insert the slot of the cap opener through the ear (picture **L**).
- Holding the bottle with one hand, turn the cap opener towards the outside of the cap (picture **M**) in order to roll the first ear onto the cap opener, then slide the ejection device (picture **N**) to eject the scrap. Repeat these operations with the second ear.
- Finally, remove completely the remainder of the side scraps with the top of the tool (picture **O**).

6- CRIMPING TOOL CHARACTERISTICS:

See technical data sheet FT 9641 and list of features RC 1279

IMPORTANT NOTE: The technical characteristics of this crimping tool and the user's guide are only provided as general information and can be updated without prior notice. TOURNAIRE S.A. will not be liable for any incorrect use of its containers and accessories or for any wrong understanding of this notice. In any case, the final user, the filling company and/or the marketing company remain the sole responsible to validate its filling and closing process, to operate according to Good Manufacturing Practices and to carry out appropriate controls of the finished product.

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