
USER'S GUIDE

T.802 closing system

IMPORTANT: This crimping tool is pre-set and tested in our premises before despatch. Any major adjustment or dismantling should be made at TOURNAIRE. In case of any doubt, do not hesitate to contact us or send us the tool for adjustment. During the 2 year warranty period, TOURNAIRE can ensure the repair or adjustment of the tool, provided it has not been submitted to abnormal mechanical or chemical constraints nor used improperly. We recommend for instance not to sterilize/ depyrogenize with oxidant substances (ethylene oxide, hydrogen peroxide,...) and rather use autoclave or Dry Heat Sterilization for this purpose.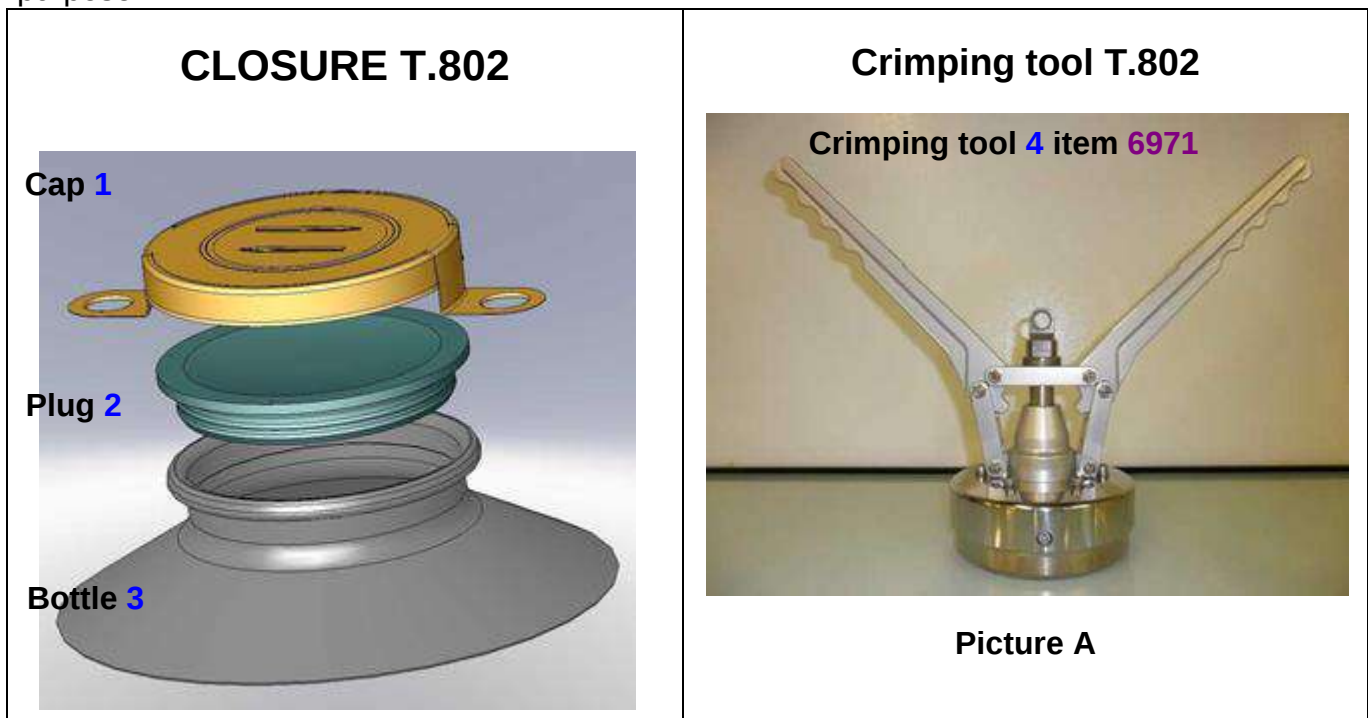

1 - PREPARATION:

1.1 Check the crimping tool before use:

Before crimping, make sure that the crimping tool 4 alone operates properly by pushing downwards the levers: you should not feel any rubbing, unusual resistance nor abnormal noise. Check also internal surface of the tool and rotating parts: they should be free from foreign objects that could impair the crimping.

In case of malfunction, please check that proper maintenance has been carried out – see **MAINTENANCE FORM NOTI-0032** or contact **TOURNAIRE S.A.** for assistance.

1.2 Preparation of bottle:

Precaution: Avoid direct contact with accessories and protect yourself with suitable gloves.

- Place the container on a flat surface at a convenient height for the operator (neck of bottle ideally at hips height).

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
2 - INSERTION OF PLUG 2:

Caution: Aluminium is much harder than the chlorobutyl stopper. Therefore during manual application, avoid slanted or folded plugs and always push gently the stopper against the aluminium neck in order to prevent risk of abrasion. Use of insertion tool is therefore highly recommended – see 2.2.

2.1 Manual insertion:

Place gently the plug **2** flat on the container neck – see picture **B Starting Phase**.

Picture B

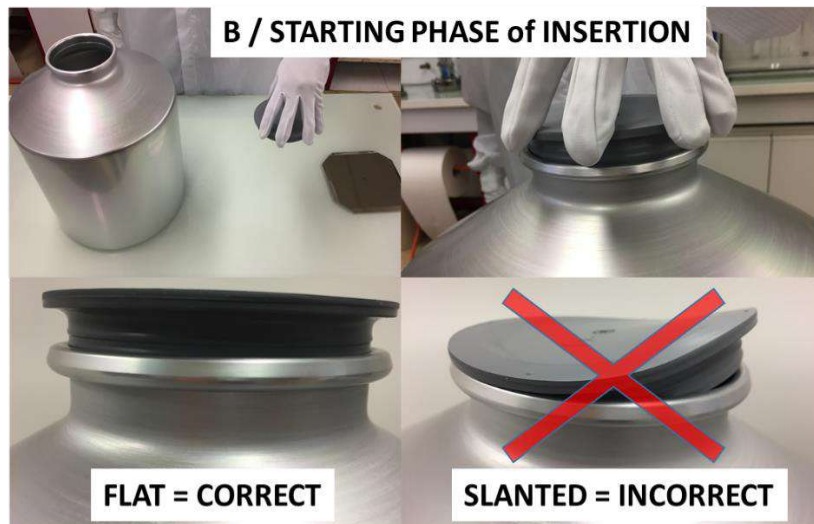Place both hands over the plug **2** and press regularly downwards – see picture **C**. Finish insertion by gently pressing all around the edge - see picture **D**. The plug must be in full contact all around the container rolled-on neck.


Troubleshooting: In case of slanted plug or rubbing or too much effort generated, check that plug is not damaged and restart insertion or discard both plug and container if plug is damaged.

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
2.2 Insertion of Plug 2 with Plug InsertionTool item 7319 (recommended way):

Important Notice: We strongly recommend using Plug Insertion Tool Ref. TOURNAIRE 7319. It indeed allows to gently bend the edge of the plug towards the inside, thus avoids the risk of abrasion, reduces the effort and simultaneously prevents potential contamination of the stopper from the gloves.


- Picking up the plug: Place the tool 5 over the plug 2. Lower the levers – see pictures E and F: plug is being sucked by the tool and can be lifted and brought to the container neck maintaining the levers horizontal – see picture G.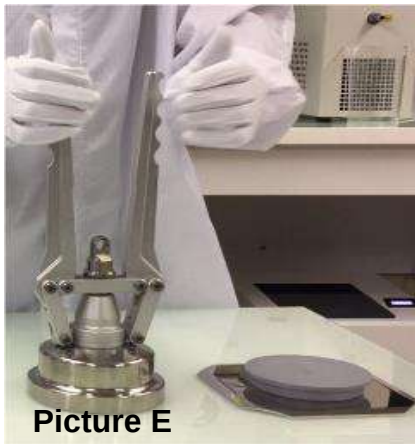

Picture E

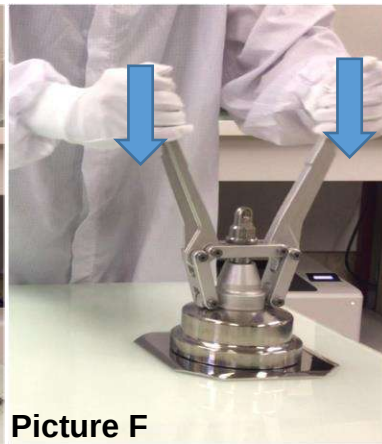Picture F

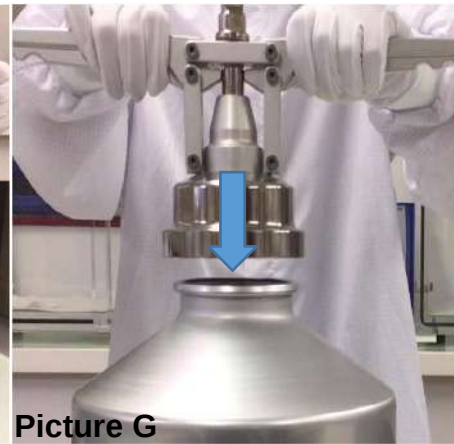Picture G

Insertion tool 5
item 7319

- Insertion of the plug: Place the tool 5 completely over the neck. Push Downwards to allow full contact of plug 2 against the neck- see picture H. Then lift the levers to release and take the tool aside- see picture I: plug is well seated inside the neck.

Picture H


Picture I

- Control of insertion: Check that the plug 2 is everywhere in full contact all around the container rolled-on neck.

3 - PLACING ALUMINIUM CAP before SEALING:


Place the aluminium cap **1** on the plug **2** on the top of it (pictures **J**, **K**, **L**), push firmly on the top, make sure that the top of cap covers entirely the plug and is in contact with the plug everywhere.

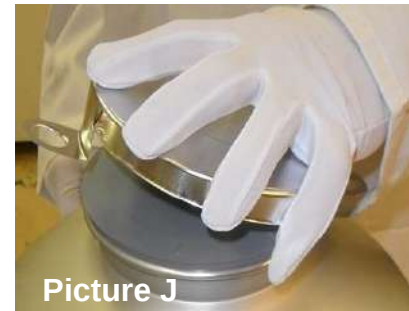Picture J

Picture K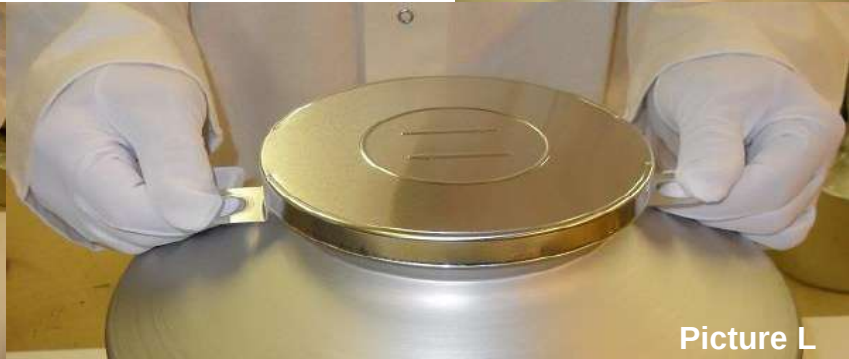

Picture L

4 - CRIMPING OPERATION:


- Position the crimping tool **4** on the assembled closure and bottle. Press firmly downwards so that the tool is strictly horizontal and fully in contact with the aluminium cap (picture **M**).
- Simultaneously push the levers downwards (picture **N**) until you reach the horizontal blocking position and levers are stopped (picture **O**). The cap is crimped.
- Lift the levers.
- Remove the crimping tool.

Picture M

Picture N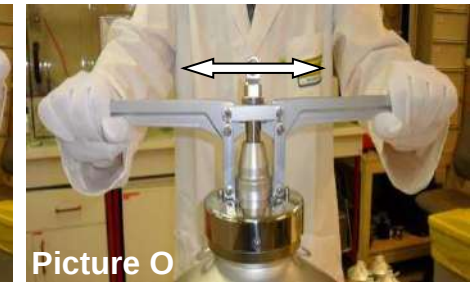

Picture O

5 - QUALITY CONTROL of the CRIMPING:

The aluminium cap **1** must be entirely rolled back on the container's neck all around the neck circumference, some little folds on the edge of the cap don't reduce the performance of the sealing (picture **P**). It is normal to see little folds on the edge of the cap.

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6 - TROUBLE SHOOTING:

In case a defective crimping is noted, first check that plug, cap and tool are properly adjusted – see § 1-1 and 1-2. If cap is blocked or deformed, take it off and use a new one then retry.

1/ too much effort is required for crimping:

- Check the crimping tool is free from any foreign object or dirt. Clean with a tissue or cloth, dipped into alcohol for instance and remove the dirt /object.
- Check that all preparation steps (§1.2) have been thoroughly respected, particularly the full horizontal insertion of the plug and cap, everywhere in contact.
- Repeat crimping operations (see - §2 -) with a new aluminium cap **1**
- Check that the crimping tool has been well maintained (see especially NOTI-0032 §4)



2/ crimping still not conform (Cf. § 3):

In case of malfunction of the crimping tool, doubt or incorrect crimping, please check that proper **MAINTENANCE FORM-NOTI-0032** has been carried out

In case of doubt, do not hesitate to contact TOURNAIRE.

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- **7 - CAP OPENING INSTRUCTION:** If not easily reachable, lift one of the cap ears upward with the edge of the cap opener **6** and insert the slot of the cap opener through the ear (picture **Q**).
- Holding the bottle with one hand, turn the cap opener towards the outside of the cap (picture **R**) in order to roll the first ear onto the cap opener, then slide the ejection device (picture **S**) to eject the scrap. Repeat these operations with the second ear.
- Finally, remove completely the remainder of the side scraps with the top of the tool (picture **T**).

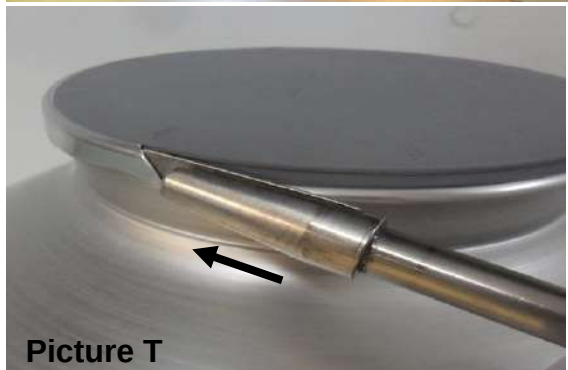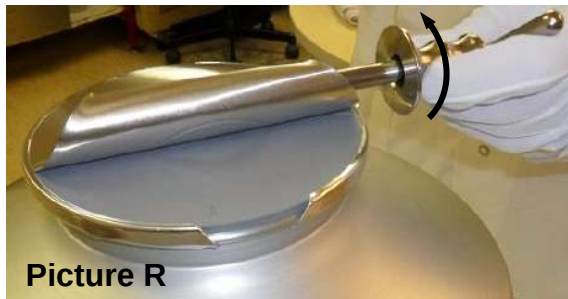IMPORTANT NOTE: The technical characteristics of this tools and the user's guide are only provided as general information and can be updated without prior notice. TOURNAIRE S.A. will not be liable for any incorrect use of its containers and accessories or for any wrong understanding of this notice. In any case, the final user, the filling company and/or the marketing company remain the sole responsible to validate its filling and closing process, to operate according to Good Manufacturing Practices and to carry out appropriate controls of the finished product.

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